

Gas Chromatograph

chrozen GC

This Value Is Not Negotiable




This Is Enough!

chro~~zen~~ GC

With accuracy and precision for each analysis, we will make a way.

YOUNG IN Chromass producing reliable and superior chromatographs for over 30 years has proudly launched our 6th generation of GC, the new ChroZen GC, by our commitment to maximize customer satisfaction with our instruments. ChroZen GC equipped with the state-of-the-art technology in both hardware and software enhances the user convenience and the data reliability while delivering the exceptional analytical solutions.

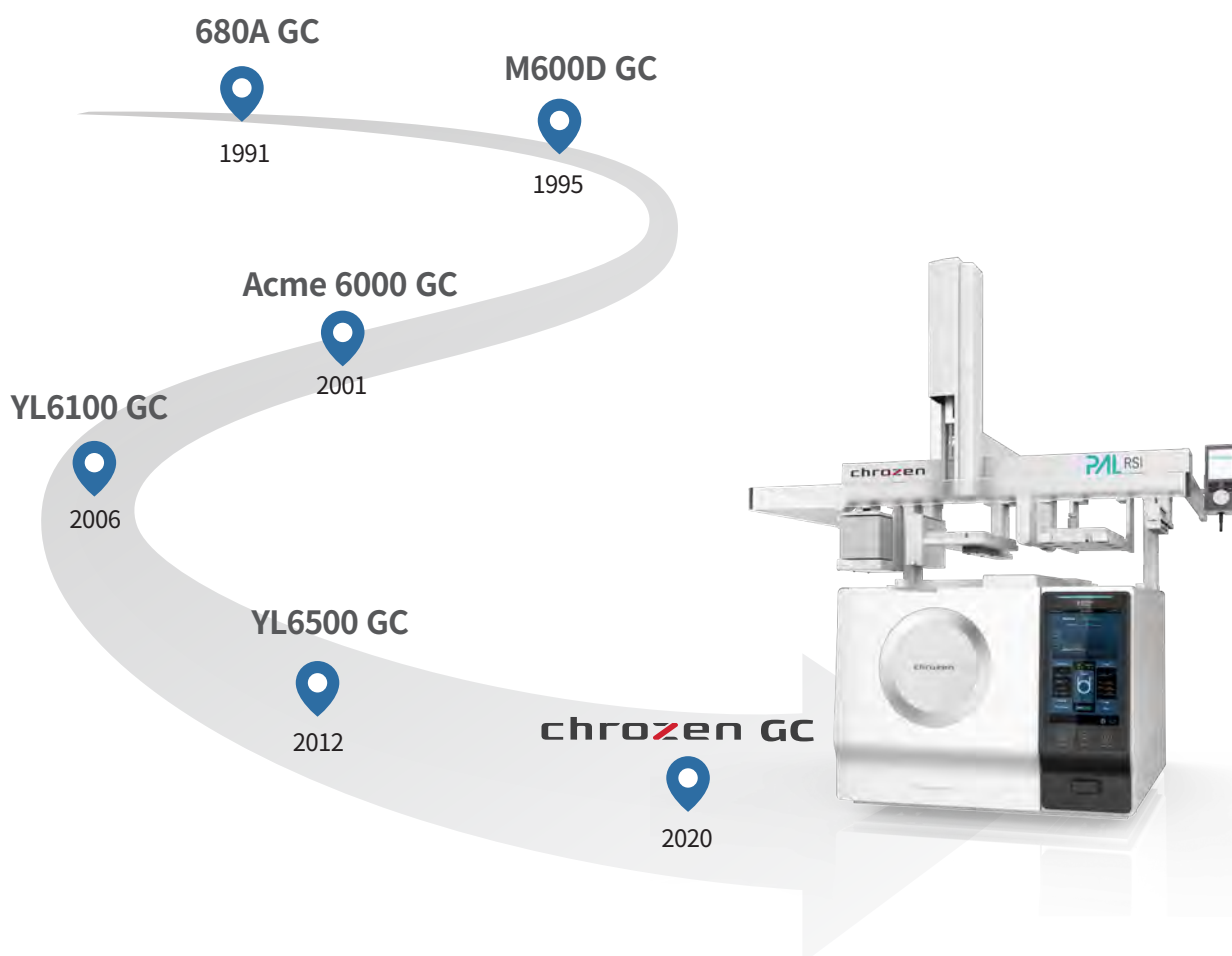

chro~~zen~~ is a brand name of YOUNG IN Chromass products that stands for the combination of **chromatograph** and **zenith**.

chro~~zen~~ GC at a Glance

The brand new ChroZen GC with its fancy design and intuitive LCD display features substantially enhanced sensitivity and reproducibility by its **powerful UPC (Ultimate Pneumatic Control)** to enable more reliable and accurate data.

More Than Smart Things

Intuitive touchpad LCD display (10.1") monitoring equipped modules (inlets, detectors, oven) at a glance and chromatogram as well as temperature programming in real time

More Than Innovation

- Substantially **enhanced sensitivity** by the efficient EMI (Electromagnetic Interference) shielding and other robust electronic parts
- **Pulsed pressure mode** (Split/Splitless): Instantly increased forefront pressure at injection minimizes the sample dispersion to provide significant sensitivity
- **Various detectors with superior sensitivity available**

More Than Reliability

- Powerfully designed UPCs (Ultimate Pneumatic Control) to optimize carrier gas control to provide exceptional accuracy and precision (**RT Repeatability SD < 0.0008 min**)
- 15 heated zones to achieve optimized thermal stability for each module (**Temperature Stability < $\pm 0.01^{\circ}\text{C}$**)
- Ultra-high speed motor to speed up the oven cooling down time for improved productivity (**$450^{\circ}\text{C} \rightarrow 50^{\circ}\text{C}$: < 3.8 min**)

More Than Solution

- Various dedicated analyzers fully configured (Up to 9 UPCs) for food, pharmaceutical, petrochemical and environmental application field that are supporting numerous standard methods such as ASTM D1945, ASTM D2163, ASTM D4059, ASTM D4815/5580, ASTM D5623, ASTM D6730, ASTM D7823 and ASTM D8071
- Extended inlet pressure range (~150 psi)
- Safe use of Hydrogen as carrier gas for superior productivity and sensitivity by H₂ leak detector (Option)
- SmartFlow Kit for backflush, 2D GC or heart-cut application to optimize analytical efficiency in multiple way



Accuracy and Precision is what ChroZen GC values most.


ChroZen GC with powerful and stable UPCs (Ultimate Pneumatic Control) ensures data reliability and facilitates the most remarkable repeatability in both RT (Retention Time) and area that considered as one of significant factors for estimating how good the GC performs.

RT(min)	C12	C14	C16
1	2.6125000	3.7200000	4.8750000
2	2.6125000	3.7200000	4.8750000
3	2.6125000	3.7208333	4.8758333
4	2.6125000	3.7208333	4.8750000
5	2.6125000	3.7200000	4.8750000
6	2.6125000	3.7208333	4.8741667
7	2.6125000	3.7200000	4.8750000
8	2.6133333	3.7208333	4.8758333
9	2.6133333	3.7208333	4.8750000
10	2.6133333	3.7216667	4.8750000
11	2.6133333	3.7200000	4.8750000
12	2.6133333	3.7208333	4.8750000
13	2.6133333	3.7208333	4.8758333
14	2.6125000	3.7208333	4.8750000
15	2.6133333	3.7208333	4.8758333
SD(min)	0.000416	0.000478	0.000451

RT Repeatability SD < 0.0008 min


Area(pA.s)	C12	C14	C16
1	0.3732875	0.3704902	0.3615789
2	0.3768552	0.3718810	0.3651483
3	0.3745179	0.3717736	0.3644548
4	0.3759930	0.3721659	0.3660563
5	0.3745244	0.3715350	0.3663568
6	0.3743780	0.3714631	0.3644980
7	0.3760052	0.3735342	0.3655185
8	0.3752988	0.3718196	0.3654820
9	0.3746218	0.3712330	0.3646219
10	0.3753269	0.3724284	0.3642208
11	0.3759125	0.3734419	0.3666292
12	0.3747429	0.3713757	0.3645661
13	0.3733229	0.3689732	0.3628517
14	0.3762151	0.3725241	0.3656627
15	0.3767046	0.3730961	0.3653833
%RSD	0.285095	0.301063	0.348607

Area Repeatability RSD < 0.5%

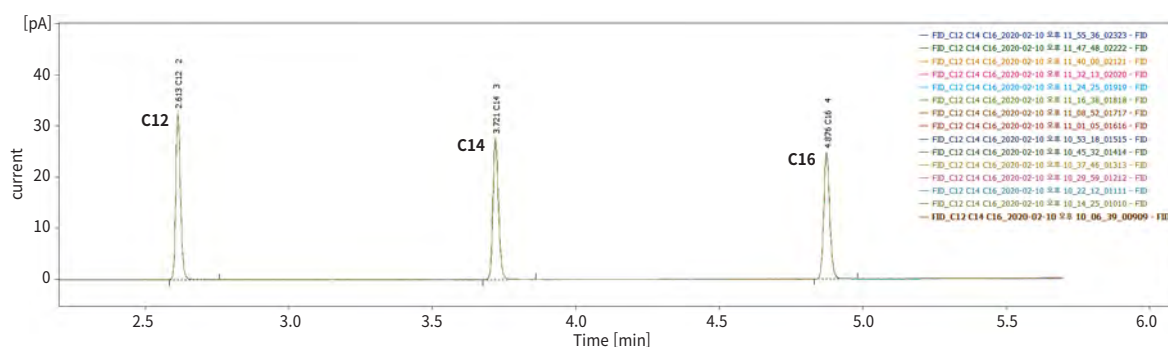


< RT Repeatability and Area Repeatability Data in 15 Sequence Injections >

Pressure Stability with Temperature Change $\leq \pm 0.002$ psi

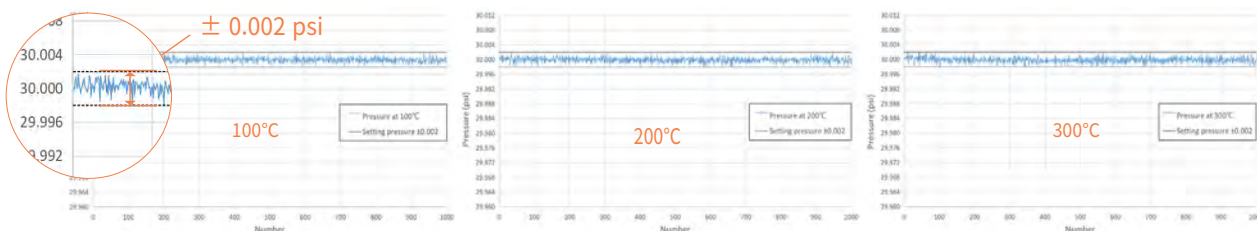

Flow Stability with Temperature Change $\leq \pm 0.001$ mL/min


- column flow rate = 4 mL/min
- measurement time = 10 min

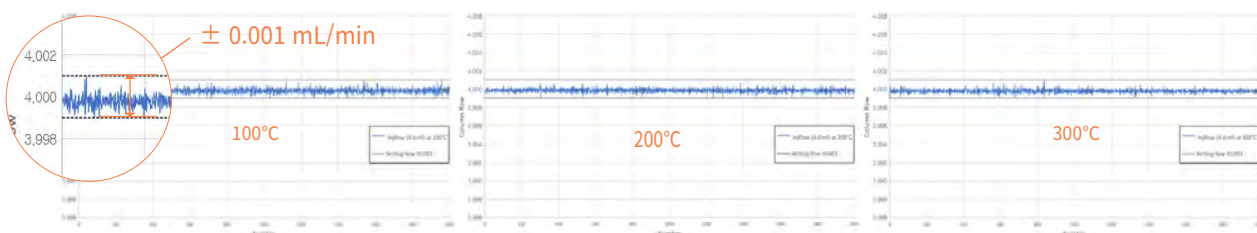

Nothing More Valuable than Your Time, ChroZen GC

Faster, The faster it cools down, the more time you save.


Ultra-high speed motor enables to speed up the oven cooling down time to reach within 3.8 minutes from 450°C to 50°C, which leads users to reduce the analysis time to improve productivity.

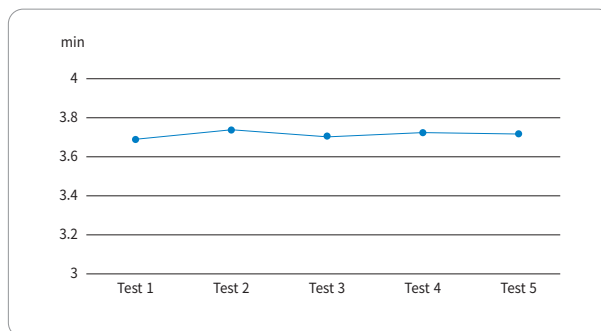

<450°C→50°C Cooling Time Test>

Much Faster, 50~70% decrease in the average analysis time with H₂ as a carrier gas!


The Van Deemter equation describes the column efficiency varying with mobile phase velocity and hydrogen has much greater velocity range while providing the desired separation and optimizing overall productivity. Thus, using a hydrogen as a carrier gas can lead significant time saving and the faster linear velocity eluted components faster that ends up with more narrow peaks, higher sensitivity in other words, while maintaining peak separations. However, all of these benefits should come with safety, which can be ensured by ChroZen GC equipped with hydrogen leak detector (Option) to avoid much of the risk.

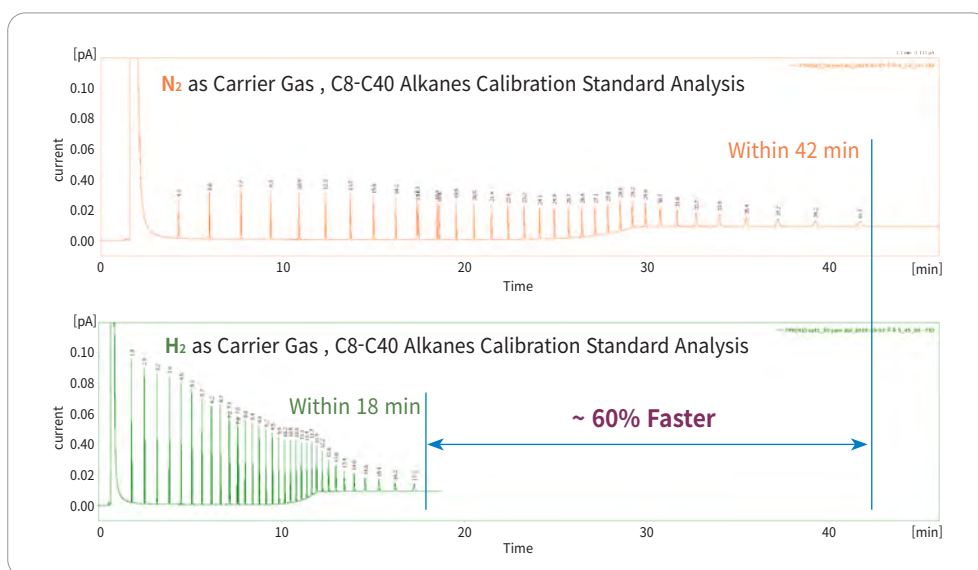

<Comparison of Hydrogen and Nitrogen as a Carrier Gas>

Much Longer, Consider target compounds only for column longevity by backflush


Backflush technique delivers the decrease in analysis time as well as the enhanced column life time by reversing the column flow after target compounds are eluted to cut the unnecessary procedure. ChroZen GC can be simply configured by adding the complete module and it gives very precise and accurate flow control for the remarkable reproducibility.

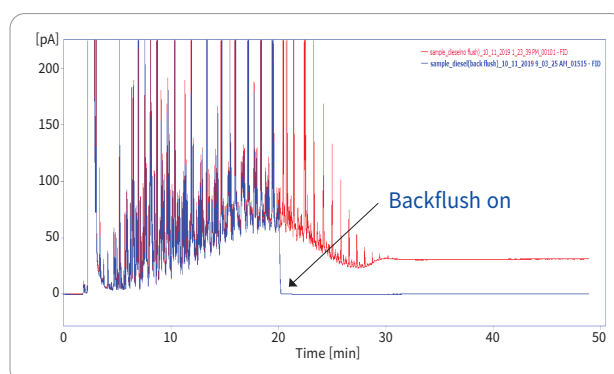

ChroZen GC Is on the Ball, Superior Sensitivity

Pulsed Pressure Mode (Available with Capillary Inlet)

In the pulsed pressure mode, ChroZen GC instantly increases forefront pressure at injection to avoid the sample dispersion and minimizes backflash exceeding the liner volume as well as carry-over to provide significant sensitivity.

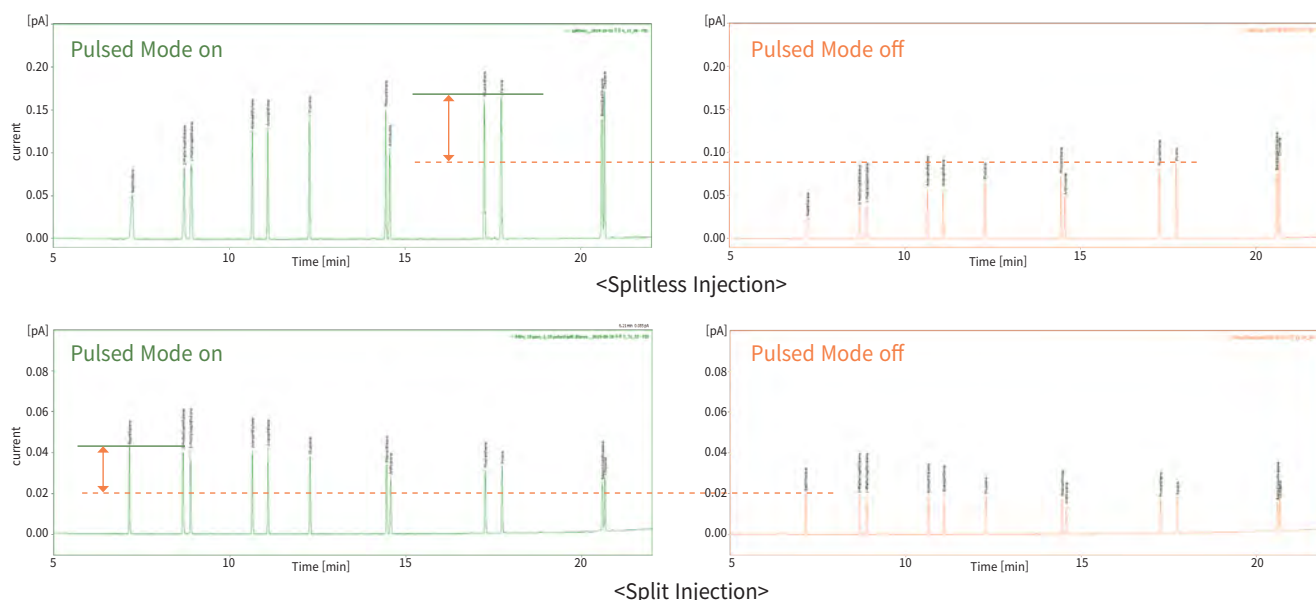


Enjoy the Variety, Remarkable Detectors

There are various detectors with superior sensitivity available for users' demands depending on their applications to make the analysis results more efficient and accurate.

- Flame Ionization Detector (FID)
- Micro-Thermal Conductivity Detector (μ TCD)
- Pulsed Discharge Detector (PDD)
- Flame Photometric Detector (FPD)
- Photoionization Detector (PID)
- Thermal Conductivity Detector (TCD)
- Nitrogen Phosphorus Detector (NPD)
- Pulsed Flame Photometric Detector (PFPD)
- Vacuum Ultraviolet Detector (VUV)

FID (Flame Ionization Detector), as one of these detectors, which is widely used for detection of hydrocarbons ensures extraordinary sensitivity with less than 1.2 pg C/s MDL (Minimum Detection Limit). After conducting thousands times of sequential tests for baseline noise and drift, our FID remains stable and reproducible data, which enables data reliability.

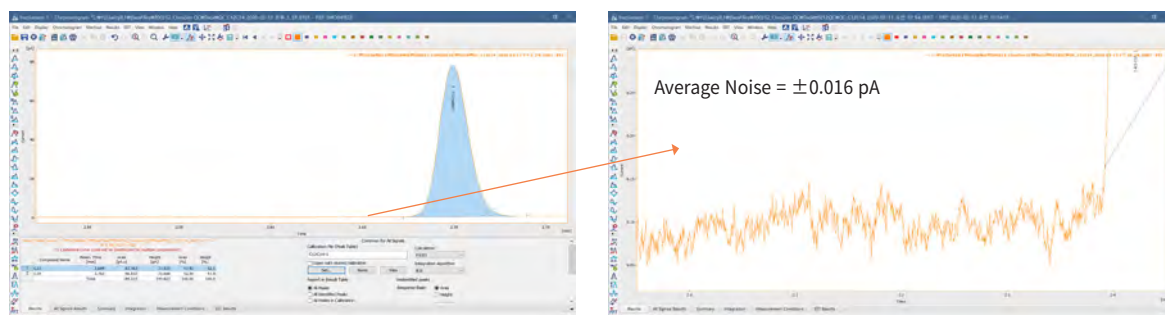

<MDL of FID in Real Data, 1.05 pg C/s>


Judge a Book by Its Cover!

Stuffy Laboratory, Get Beauty


ChroZen GC with its fancy design is equipped with the intuitive touchpad LCD display (10.1"), which enables easy control of all the installed modules (inlets, detectors and oven) and monitors chromatogram as well as temperature programming steps in real time to control whole process at a glance.



High Resolution Vivid LCD with Intuitive GUI

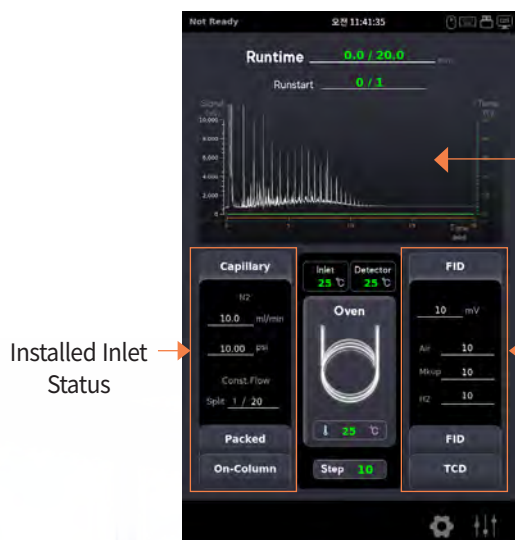

Chromatogram Monitoring in Real Time

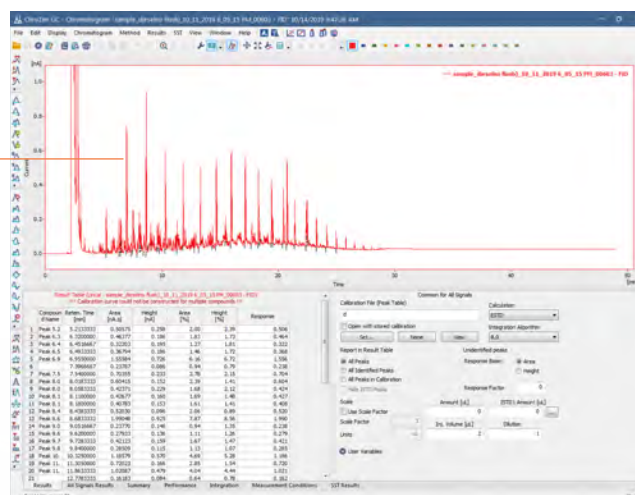

Installed Inlet Status

Installed Detector Status




Of the User, By the User, For the User

Dedicated Analyzer

Our accumulated experience in gas chromatography and devotion for customers' satisfactions led to supply one stop solution for various dedicated applications. You only need to let us know what to analyze, and then every single component that is required to analyze your sample will be configured right away.

Natural Gas Analyzer

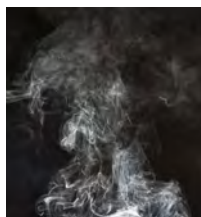

Natural Gas Analyzer with accurate valve switching function vents unnecessary components and sends significant things in time to the appropriate columns and detectors for optimized analysis of natural gas containing low level of oxygen (< 3%), isomers and compounds more than C4 which are not needed for analysis results.

* Standard Method Support: ASTM D1945, etc.

Refinery Gas Analyzer



With the availability of no. of installation (3 for each detector and inlet), Refinery Gas Analyzer can configure several columns, switching and sampling valves as well as the appropriate detectors to analyze the complex and difficult refinery gas samples.

* Standard Method Support

: ASTM D4815, D5580, D7423, D6730, D1946, etc.

Residual Solvent Analyzer



Residual Solvent Analyzer coupling with headspace autosampler can accurately and efficiently detect and quantify residual solvents found in the inks used for the printing of packaging and raw materials for food and drug products, which are known to be hazardous to human health if ingested.

Residual Pesticide Analyzer



Residual Pesticide Analyzer equipped with ECD or NPD is the optimized system to separate complex mixture of compounds containing organophosphorus and organochlorine pesticides and to run EPA methods. We also provides the analyzer equipped with GC/MS for quantitation of pesticides at trace levels.

Fatty Acid Analyzer



Fatty Acid Analyzer is suitable to analyze various fatty acids to determine the ratio of saturated fatty acids and unsaturated fatty acids in fat and provides all the solution from the preparation to the method set-up.

Phthalate Analyzer



Phthalate Analyzer excels at detection and quantitation of regulated phthalates that became concern and into force in restrictions on the use of certain substances such as children's toys, medical devices and food packaging, which are directly related to the human health.

VOC Analyzer



VOC analyzer optimized to analyze VOCs in aqueous samples that can have a bad long-term influence on human health accurately analyzes trace level of VOCs (ppt level by P&T or SPME Arrow). It equips all necessary configuration for the analysis including detailed procedures and protocols, which are in full compliance with approved U.S. EPA methods.

Biodiesel Analyzer



Biodiesel Analyzer identifies and quantifies impurities in biodiesel as an alternative fuel today and provides all solution including the sample preparation and the method set-up while achieving excellent reproducibility and resolution.

Sulfur Analyzer



Sulfur Analyzer provides increased sensitivity and selectivity with PFPD (Pulsed Flame Photometric Detector) for the analysis of sulfur compounds at trace levels as well as the analysis method achieving ASTM 6228.

Save Your Time in Laboratories

Liquid Autosampler



ChroZen PAL LSI

ChroZen PAL LSI

ChroZen PAL LSI (Liquid Sample Injection) System can reach any three-dimensional position within its working space and is ideally suited to pick and place objects like vials. It can inject from several sample containers into different inlets or transport vials from a tray to a temperature controlled modules like the peltier stack for cooling them down to 4°C.

- **Large sample capacity:** 648 position of 2 mL vials
- **Discrimination free injection:** Ultrafast injection down to 100 msec reduces needle discrimination in GC to C40/C20 > 0.98



YL3100A Autosampler

YL3000A Series Autosampler

Its sophisticated sampling techniques provide the variable needle depths to draw the sample anywhere within the vial allowing you to use various types of syringes with volume from 0.5 to 100 µL to cover the widest range of applications.

- **The self-aligning “plug and play”:** Easy positioning/Installation/re-positioning
- **Full-color touch screen interface:** Easy system accessibility and usability
- **Double wash step capability:** Pre-and post-wash for superior analytical performance (carry-over free)
- **Syringe illumination:** Easy check against air bubbles in method validation
- **Syringe ID:** Automated syringe identification and tracking of the syringe consumption based on RFID tags



YL3000A Liquid Autosampler

	YL3100A	YL3000A	ChroZen PAL LSI
Operation Type	Rack-Turret	Tray in-out and Turret	X - Y - Z Position
Sample Capacity (2 mL)	15 vials	121 vials	Max. 648 vials
Washig Vial	6 (10 mL)	6 (10 mL)	4 (10 mL) or 2 (100 mL)
User Interface	Key Pad/ Software	Touch Screen/ Software	Handheld Controller/ Software
Installation Kit	Simple bracket included already	An installation kit is requested	A mounting kit is requested
Dual Injection (In Sequence)	Available	Available	Available
Integrated Bar Code Reader	N/A	N/A	Available as Option
Syringe ID	N/A	N/A	N/A

Headspace Autosampler



YL2000H Headspace Autosampler



YL2100H Headspace Autosampler



Versa Headspace Autosampler

YL2000H Series Headspace Autosampler

The YL2000H with a heated gas-tight syringe eliminates tubing, dead volume and sample absorption for system robustness and low maintenance frequency and its compact installation on most GCs allows you to utilize headspace technology in the efficient way.


- User-friendly touch screen
- Easy installation on most GCs
- Adjustable sample volumes without loop changes

	YL2000H	YL2100H	Versa	ChroZen PAL RSI/RTC
Injection Type	Gas Tight Syringe	Gas Tight Syringe	Pressured Loop	Gas Tight Syringe
Extraction	Static HS	Static HS	Static HS	Static HS & Dynamic HS
Sample Capacity	42 vials (20 mL)	14 vials (20 mL)	20 vials (22 mL)	180 vials (10/20 mL)
Oven Position	6	1	1	6
Interface	Touch Screen /Software	Keypad /Software	Software	Hand-Held Terminal/ Software
Syringe Volume (Standard)	2.5 mL	2.5 mL	1 mL	2.5 mL
Syringe Volume (Option)	1, 5 mL	1, 5 mL	100, 250, 500 µL and 2, 3, 5 mL	1, 5mL

All in One Autosampler



YL2800T All in One Autosampler

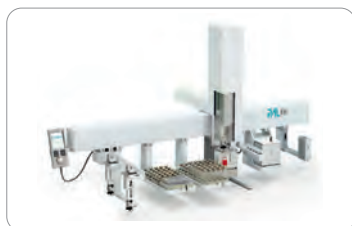

ChroZen PAL RSI/RTC

YL2800T Series Autosampler

YL2800T is a single unit combining static headspace analysis, liquid sample injection and SPME (Solid Phase Micro-Extraction) as an option.

- Sample Capacity: Headspace/SPME - 42 vials (20mL) Liquid - 121 vials (2mL)

ChroZen PAL RSI/RTC System

ChroZen PAL RSI/RTC system can be adapted or extended to provide the combined injection techniques such as static headspace injection, liquid injection, SPME (Solid Phase Micro Extraction) and ITEX (In-Tube Extraction) dynamic headspace in one instrument. ChroZen PAL RTC system provides same reliability, flexibility and productivity as the RSI system but it additionally offers the robotic tool change that switches between different tools automatically.

Applying smart syringe technology with ID chip, it automatically preset all of syringe parameters, ranges and usage tracking for users' convenience.

Sample Preparation System



Lumin Purge and Trap Concentrator

The Lumin Purge and Trap Concentrator (PTC) is a sample preparation instrument used to remove Volatile Organic Compounds (VOCs) out of aqueous and solid sample types using Helium or Nitrogen.



AQUATEk LVA Waters-only Autosampler

The AQUATEk LVA is a Purge and Trap (P&T) Autosampler that automates the sample preparation steps for the analysis of liquid samples via purge and trap. The system is capable of preparing samples such as drinking water and wastewater.



Atomx XYZ Automated VOC Sample Prep System

The Atomx XYZ is the second generation combined soil/water autosampler and purge and trap concentrator system to analyze VOCs. It provides 84-position vial tray with optional vial chiller to 4°C and the improved moisture control system reduces the amount of water transferred to the GC.



ChroZen ATD

ChroZen ATD is an automated thermal desorption system which enhances productivity by 50 sorbent tubes capacity and smart sequence overlapping function. With two-stage desorption, it also utilizes versatile thermal desorption by concentrating volatile and semi-volatile organic compounds collected on a sorbent tube into a focusing trap to enable the analysis of VOCs at trace level with the superior sensitivity.



Pyrolyzer, Pyroprobe 6000 series

Pyrolyzer coupled with ChroZen GC allows you to analyze the samples such as paint, tapes, caulk, adhesives, food packaging, rubber, plastic, papers, ink, coating and a full range of household products without extractions or derivatizations for qualitative and quantitative information.




Gas Sampling Valve/Purge Housing

In order to obtain accurate results in gas analysis, it is necessary to equip with a gas sampling valve in the front of column. The gas sampling valve enables manual or automatic gas sample injection and flow switching. Our gas sampling valve is operated by an air actuator or μ -electronic actuator which opens and closes it. The Purge housing eliminates any possible diffusion from the atmosphere into the valve, or safely vents fugitive emissions from the valve to analyze samples of O₂, N₂, etc at low concentration of a few ppms.

Powerful and Intuitive Control, YL-Clarity

The sophisticated YL-Clarity chromatography data system is easy to use and offers extensive data management plus full control of our GC products by its intuitive user interface. The software is designed for 21 CFR Part 11 Compliance and full compatibility with MS Windows OS and seamlessly handles data processing and instrument control using an ultra-reliable LAN communication.

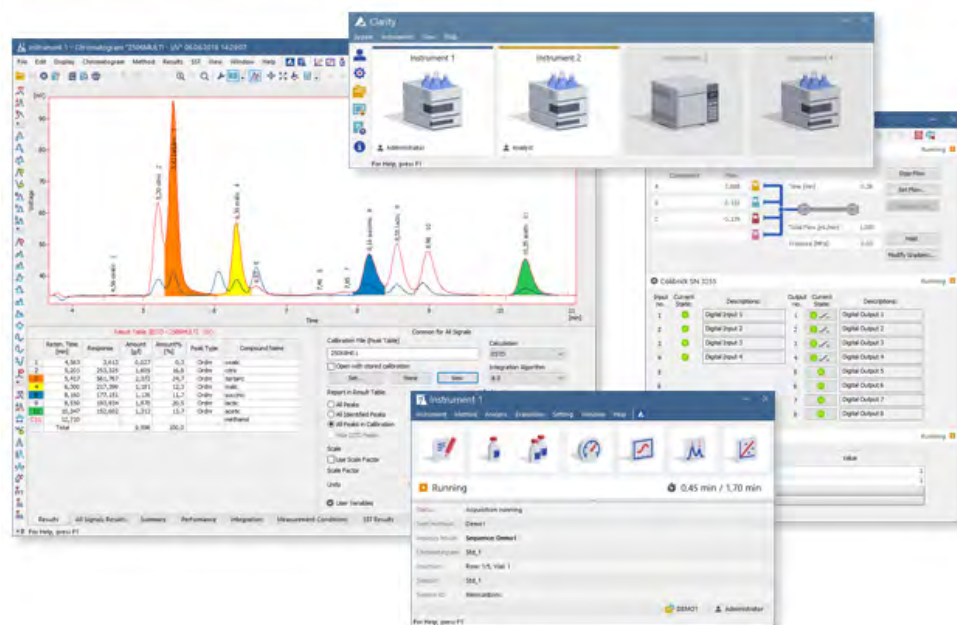

- List of Controllable Instruments: ChroZen GC, ChroZen GC/MS, YL6500 GC, YL6900 GC/MS, ChroZen UHPLC, YL9100 Plus HPLC, YL9300 HPLC, YOUNG IN Chromass autosamplers
- MS Windows Compatibility: Vista, 7, 8, 10, 11
- Optional Module
 - SST (System Suitability Test)
 - NGA (Natural Gas Analysis)
 - PDA Extension
 - GPC Extension
 - MS Control
 - DHA (Detailed Hydrocarbons Analysis)
 - Validation Kit
- 21 CFR Part 11 Compliance: Management of user accounts, audit trail and electronic signature
- Simultaneous data acquisition from up to four independent chromatographs: Each chromatograph can acquire data from up to 32 detectors.

Powerful GC/MS Library

Various GC/MS Libraries are available depending on specific applications with multiple information of each compound's spectra, chemical structure and retention index to get the accurate and reliable data.

GC - MS Data			
Library	Wiley Registry 2023	NIST 2023	WR23/NIST 2023 GC-MS Total
Spectra	873,300	394,000	1,180,000
Chemical Structures	841,100	347,100	1,148,600
Compounds	741,000	341,600	950,200



Gas Generator



Air Generator

(Zero Air-GC Series/Ultra Zero Air-GT Series)

Air Generators produce dry and hydrocarbon-free air (NO_x/SO_x free and CO_2 free with GT series) by using air from oil free compressors, thus avoid the need to use conventional bottles that are often complex to change. With this zero air, you can decrease the background noise level because it gives the baseline much better stability, considerably increasing detector sensitivity and ensuring precise analytical results.



N₂ Generator

(N₂ Tower Plus)


N₂ TOWER Plus produces high purity N₂ (up to 99.999%) for all major applications in analytical laboratories. Nitrogen generators eliminate the typical problems associated with the handling of high pressure gas tanks while reducing the cost for high purity nitrogen at the same time.

YOUNG IN Chromass

As one of the leading analytical instrument manufacturers, YOUNG IN Chromass has produced reliable and superior chromatographs over more than 30 years by dedicating our research and development to chromatographic technologies in Korea.

We provide our customers with reliable analytical instruments at reasonable prices and keep constantly collaborating with our customers and sharing with them valuable information on experiments in science, on chemical analysis and all others. In more than 50 countries, we are ready to back you up, relying on our well-trained distributors wherever and whenever new needs arise. We will provide our best support for all of our customers to select optimal instruments with a right solution to application and by expanding our valuable research and experiments.

Our Mission Is to Provide Entire Solutions for Chromatography

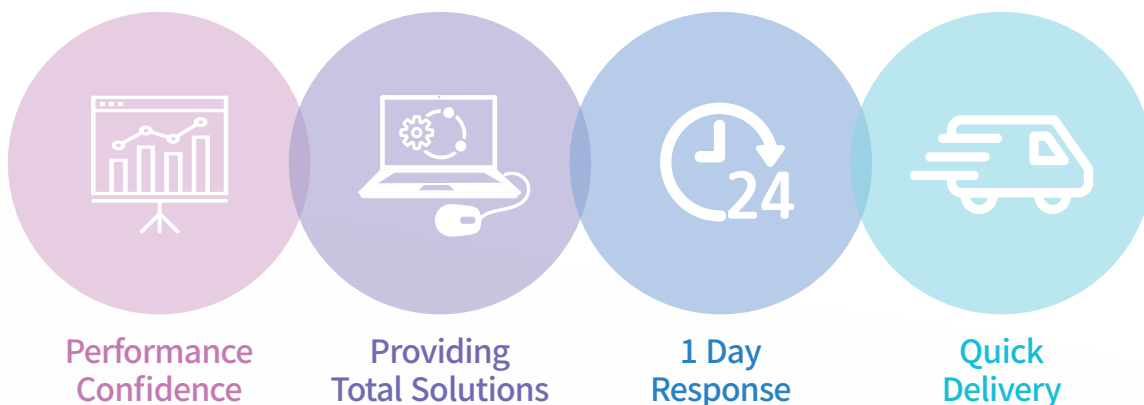

Customer Support





60, Anyangcheondong-ro, Dongan-gu, Anyang-si,
Gyeonggi-do, 14042, Korea
TEL : +82-31-428-8700 / FAX : +82-31-428-8787
E-mail : export@youngincm.com
Homepage : www.youngincm.com



Young In Chromass products are endorsed by
Korean PPS(Public Procurement Service) in
recognition of their excellent technologies and
the product quality.